


New project: semi manual bag filler BT1

für Promens Intracon and Intracon single neck infusion bags

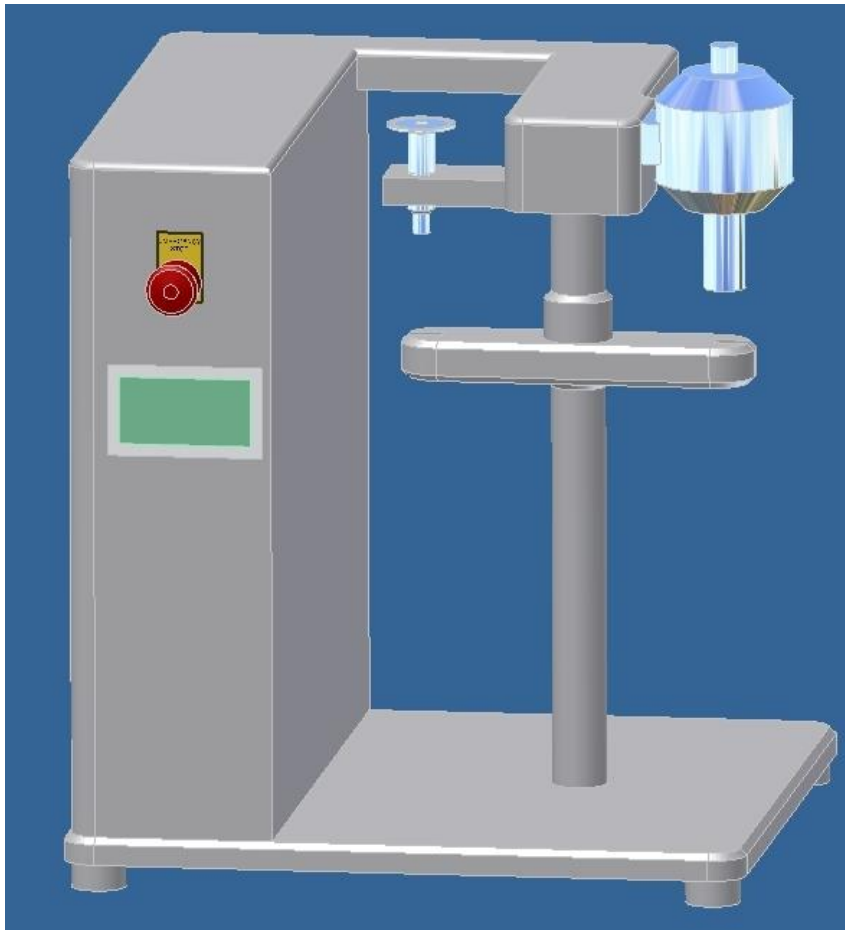

Bag types:	Intracon and Intracon single neck up to 1000 ml
Vial types:	10-100 ml
Bottle types:	100-1000 ml (with extra change parts)
Stopper and cap size:	20 mm (Standard), 32 mm (with change parts)



Technical data:

Dosing system:	massflow-meter or peristaltic dosing pump
Standard equipment:	basic unit with swivel arm filling station crimping station
Product contacting parts:	stainless steel 306L
Filling hoses:	silicone or EPDM
Sealing rings:	EPDM
Base frame:	stainless steel 304
Connecting parts:	stainless steel 304 and anodized aluminium
Interface:	Monochrome touch screen

Mode of operation:

- manual in-feed of empty bag into the swivel arm (right hand side)
- manual removal of protection cap
- after pressing the start button, the swivel arm turns by 90 degrees and the bag will be filled
- manual supply of stopper and alu cap as well as of next bag (without protection cap)
- after pressing the start button again, the swivel arm turns by 90 degrees and the bag will be sealed at the same time the next bag is filled